



Quick User guide

- Fridge-tag® 2
- Fridge-tag® 2 L

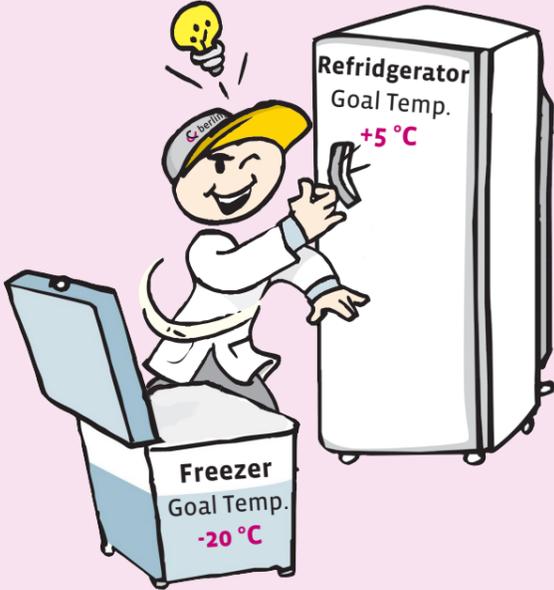

1 Reading the display during measurement

OK display

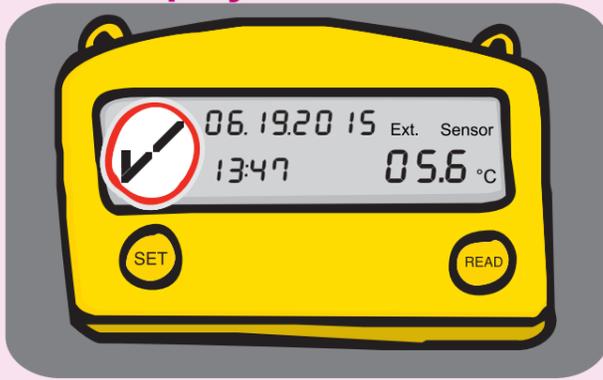

The following information is visible on screen:

- temperatures are OK
- Current date: 06.19.2015
- Current time in military format: 13:47
- Current temperature: +5.6 °C

ALARM display




If an alarm occurs the following additional symbols appear on screen:

- **ALARM** symbols, temperature alert
- = Max higher (HI) temperature ALARM
- = MIN lower (LO) temperature ALARM
- = ALARM sensor disconnected for +15 minutes

The arrow appears under today and / or the day an ALARM occurs

2 Read history (record daily checks (2x) in memory and clear an X if alarmed)

Recording daily checks

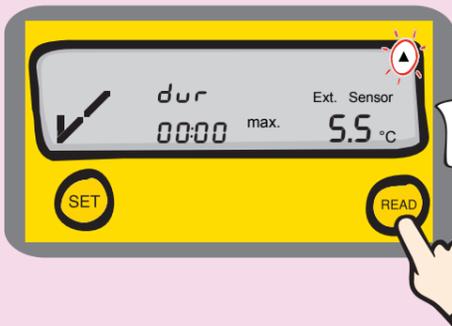


Press **READ** 3 times

- Press the READ button 3 times to record your check
- **1st** press: see UP arrow blink = HI temp. so far today
 - **2nd** press: see DOWN arrow blink = LO temp. so far today
 - **3rd** press: records in memory the time of your manual check

- A **Fridge-tag® 2** records an AM or PM for morning or afternoon checks
- A **Fridge-tag® 2 L** records the actual time it is checked in the morning and in the afternoon.

X Clearing the X / Example of a LO ALARM

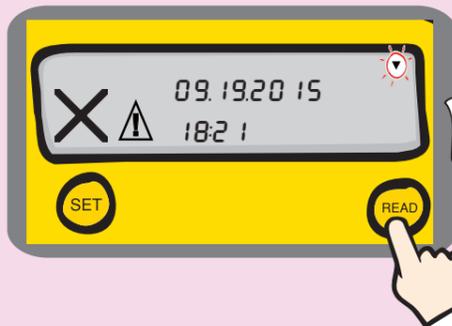

1st Press

- Press the READ button 4 or 5 times to clear the X and record your check.
- There will be an UP or DOWN arrow under today in the upper right of the screen indicating a HI or LO ALARM

If both arrows are present then it is typically an ALARM for the external sensor being disconnected for more than 15 minutes. Double arrows need 5 presses of READ to clear the X.



2nd press



3rd press




4th press

- 1st** press: see UP arrow blink = HI temp. so far today
- 2nd** press: see DOWN Arrow blink = date & time of the alarm
- 3rd** press: see DOWN arrow blink & dur = duration of ALARM and the lowest temperature during that alarm
- 4th** press: clears the X and records in memory

3 Downloading the report

STEP 1

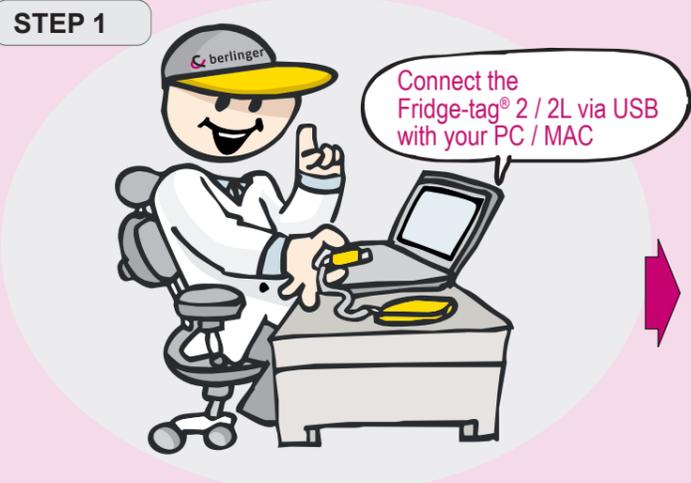

Connect the Fridge-tag® 2 / 2L via USB with your PC / MAC

Download the report from the Fridge-tag® device by connecting it to any PC / MAC.

- **Fridge-tag® 2** = 60 day report
- **Fridge-tag® 2 L** = 56 day report (with 8 weekly graphs)

STEP 2



Click on the Fridge-tag file to open it

STEP 3



Double click on the PDF file to open the report. **Note:** store reports for 3 years to be VFC compliant

4 Support



To change date / time or Celsius to Fahrenheit, view videos at: www.berlinger.com/videos